


Display & buttons

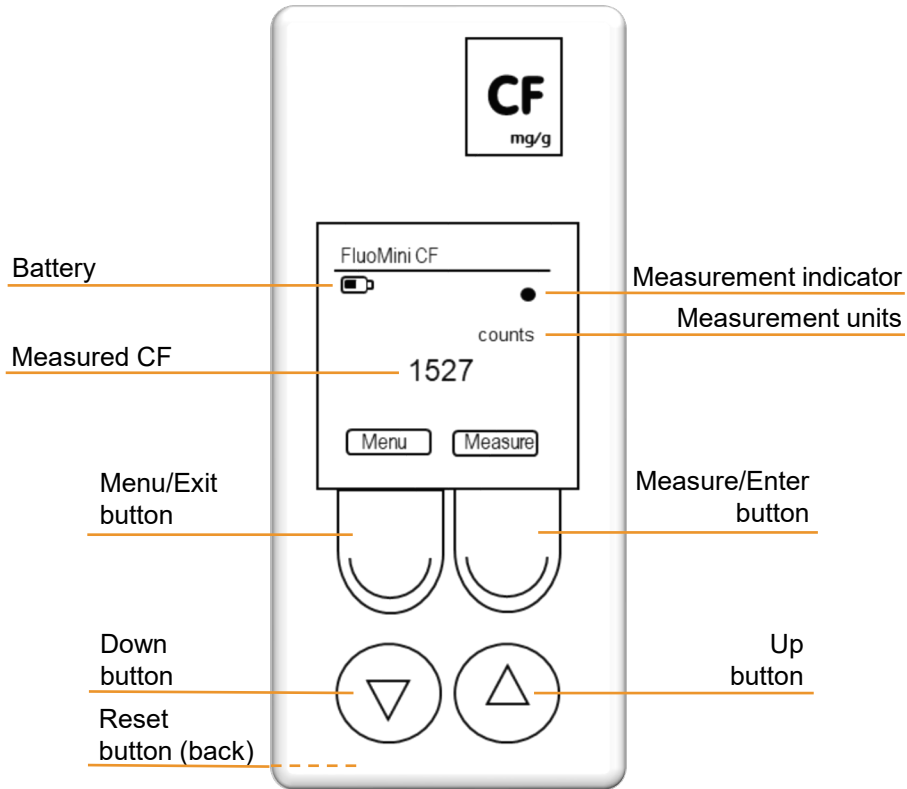

Menu/Exit button



Enter & exit main menu

Measure/Enter button



End sleep mode
Single measurement > press briefly
Start/stop continuous measurement > press 2 sec

Calibration

Apply reference sample

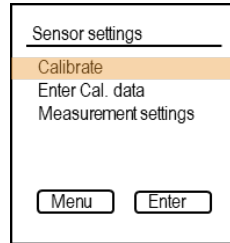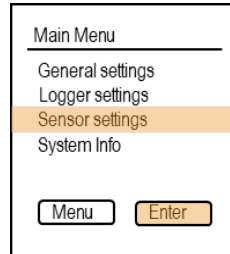




Choose between 1- or 2-point calibration

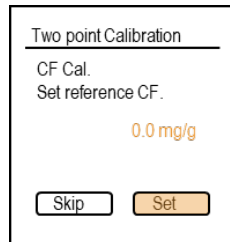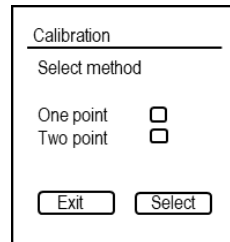




* Wait for a stable measurement

Log function

Set interval time

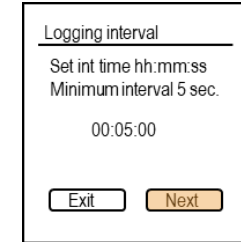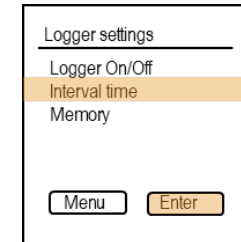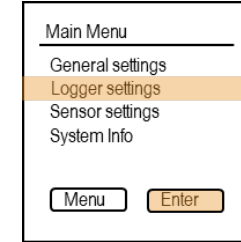



Start logging

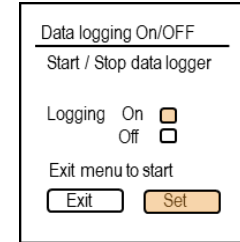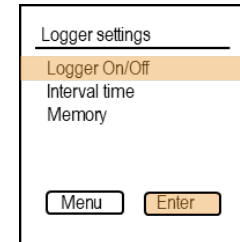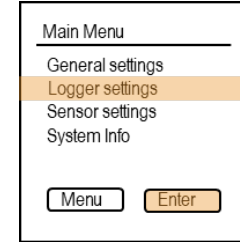

CAUTION!

When the probe touches the surface of the sample, it might get contaminated resulting in false measurements. Ensure to keep the tip of the probe clean. For further information read the manuals on our website.

Remove reference sample

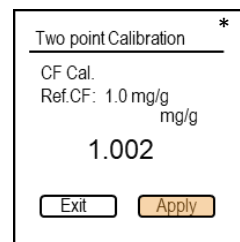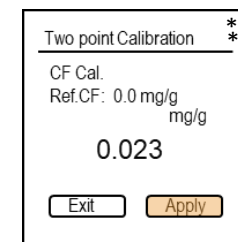

** Only for 2-point calibration